


Securitas Environmental Monitoring Custom View and Preferences Setup Tip Sheet

How to set up a custom dashboard view for Environmental Monitoring in MobileView Application:

1. In MobileView web application version 5.6 and above, we have the ability to create custom views for Environmental Monitoring. First Login to application using the web link below with your NetID Credentials:

<https://freezermonitor.its.yale.edu>


2. From the Environmental Monitoring Application Home Screen, click the three bars in the upper left to reach the Configure Views screen.

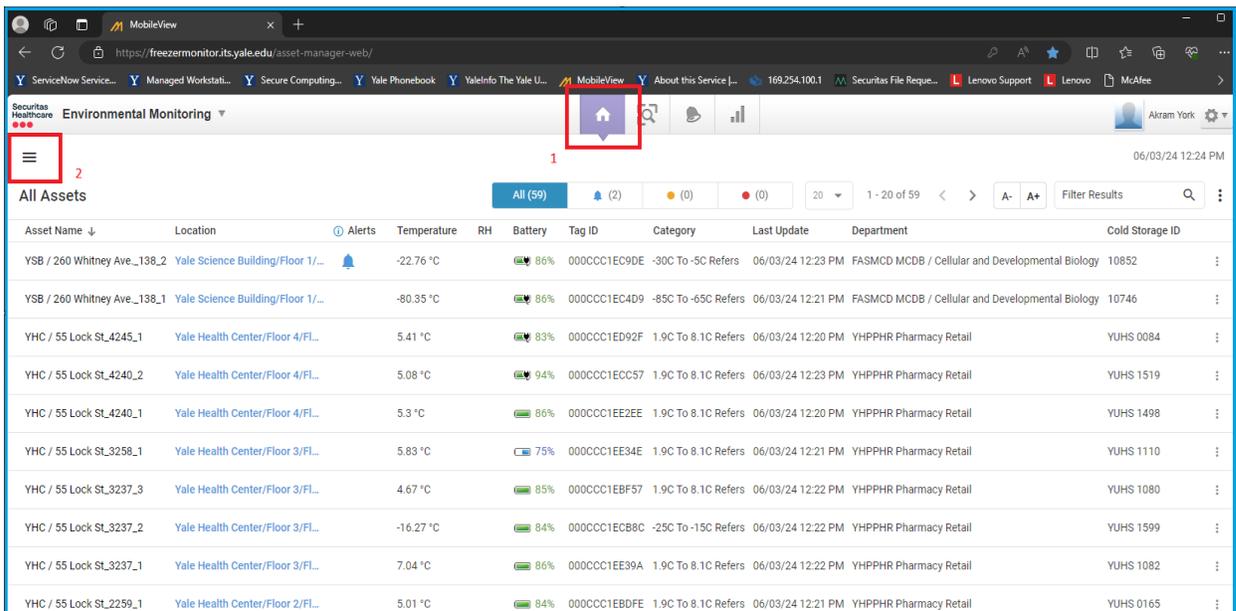

The screenshot displays the MobileView web application interface for Environmental Monitoring. The top navigation bar includes a home button (house icon) and a menu icon (three horizontal bars). The main content area shows a table of assets with the following columns: Asset Name, Location, Alerts, Temperature, RH, Battery, Tag ID, Category, Last Update, Department, and Cold Storage ID. The table lists 10 assets, each with its location, temperature, relative humidity, battery level, tag ID, category, last update time, department, and cold storage ID.


Asset Name	Location	Alerts	Temperature	RH	Battery	Tag ID	Category	Last Update	Department	Cold Storage ID
YSB / 260 Whitney Ave...138_2	Yale Science Building/Floor 1/...		-22.76 °C		86%	000CCC1EC9DE	-30C To -5C Refers	06/03/24 12:23 PM	FASMCD MCDB / Cellular and Developmental Biology	10852
YSB / 260 Whitney Ave...138_1	Yale Science Building/Floor 1/...		-80.35 °C		86%	000CCC1EC4D9	-85C To -65C Refers	06/03/24 12:21 PM	FASMCD MCDB / Cellular and Developmental Biology	10746
YHC / 55 Lock St_4245_1	Yale Health Center/Floor 4/FL...		5.41 °C		83%	000CCC1ED92F	1.9C To 8.1C Refers	06/03/24 12:20 PM	YHPPHR Pharmacy Retail	YUHS 0084
YHC / 55 Lock St_4240_2	Yale Health Center/Floor 4/FL...		5.08 °C		94%	000CCC1ECC57	1.9C To 8.1C Refers	06/03/24 12:23 PM	YHPPHR Pharmacy Retail	YUHS 1519
YHC / 55 Lock St_4240_1	Yale Health Center/Floor 4/FL...		5.3 °C		86%	000CCC1EE2EE	1.9C To 8.1C Refers	06/03/24 12:20 PM	YHPPHR Pharmacy Retail	YUHS 1498
YHC / 55 Lock St_3258_1	Yale Health Center/Floor 3/FL...		5.83 °C		75%	000CCC1EE34E	1.9C To 8.1C Refers	06/03/24 12:21 PM	YHPPHR Pharmacy Retail	YUHS 1110
YHC / 55 Lock St_3237_3	Yale Health Center/Floor 3/FL...		4.67 °C		85%	000CCC1EBF57	1.9C To 8.1C Refers	06/03/24 12:22 PM	YHPPHR Pharmacy Retail	YUHS 1080
YHC / 55 Lock St_3237_2	Yale Health Center/Floor 3/FL...		-16.27 °C		84%	000CCC1ECB8C	-25C To -15C Refers	06/03/24 12:22 PM	YHPPHR Pharmacy Retail	YUHS 1599
YHC / 55 Lock St_3237_1	Yale Health Center/Floor 3/FL...		7.04 °C		86%	000CCC1EB39A	1.9C To 8.1C Refers	06/03/24 12:22 PM	YHPPHR Pharmacy Retail	YUHS 1082
YHC / 55 Lock St_2259_1	Yale Health Center/Floor 2/FL...		5.01 °C		84%	000CCC1EBDFE	1.9C To 8.1C Refers	06/03/24 12:21 PM	YHPPHR Pharmacy Retail	YUHS 0165

3. From the Configure Views sidebar click the “+CREATE VIEW” Link.

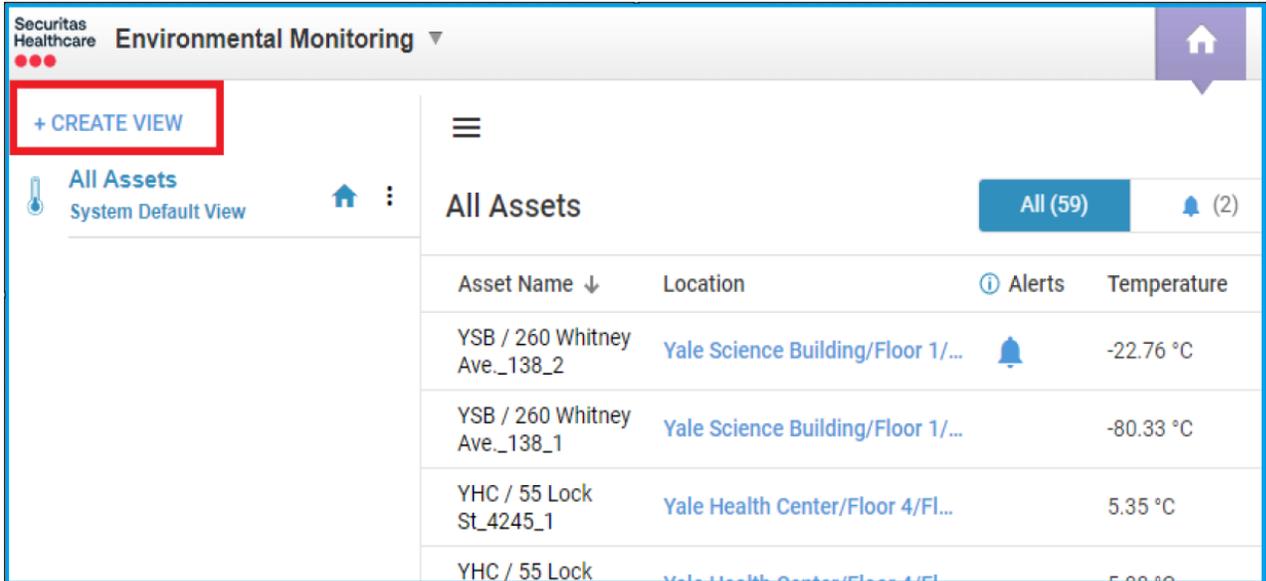

4. The Create View menu will appear on the right-hand side of the screen. Give the view a name and optional description. In this example we will create a view for the Test Department.

In the Data Filter Section, select Departments. Then scroll down and search for the Test Department and select it.

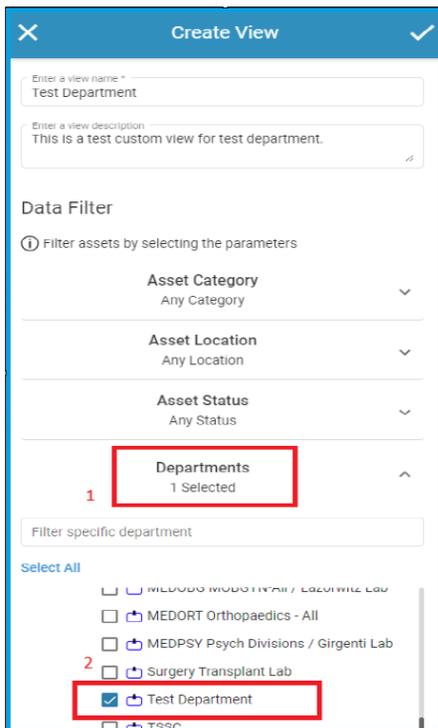


5. Scroll down to the Edit Columns Section and select the desired columns. You may also reorder the selected columns by clicking the Order Columns button. Selected Columns shown below are recommended for Environmental Monitoring Users.
6. Click the Check sign in the upper right to save the configuration.

The screenshot shows the 'Create View' dialog with the 'Edit Columns' section. The 'Select Available Columns' tab is active. A list of columns is shown with checkboxes. The following columns are checked: Asset Name, Temperature, Alerts, Department, Battery, Location, RH, Tag ID, and Last Update. Other columns like Custom Sensor, Business Status, Last BD Session, BD Status, Voltage, and various Alert_Notify and Cold Storage ID options are unchecked.

The screenshot shows the 'Create View' dialog with the 'Edit Columns' section. The 'Order Columns' tab is active. A list of columns is shown in a vertical stack, each with a three-dot menu icon to its right. The columns are ordered as follows: Asset Name, Location, Temperature, RH, Alerts, Battery, Tag ID, Department, and Last Update. Below this list is the 'General Settings' section, which includes a label 'Select location path to display:' and a row of four buttons: {BUILDING} ::, {FLOOR} ::, {AREA} ::, and {ZONE_STR} ::.

7. To make this your default view, Click the three bars in the top left of the screen again to access the View sidebar. Click the 3 dots next to your View and click "Set as Default".

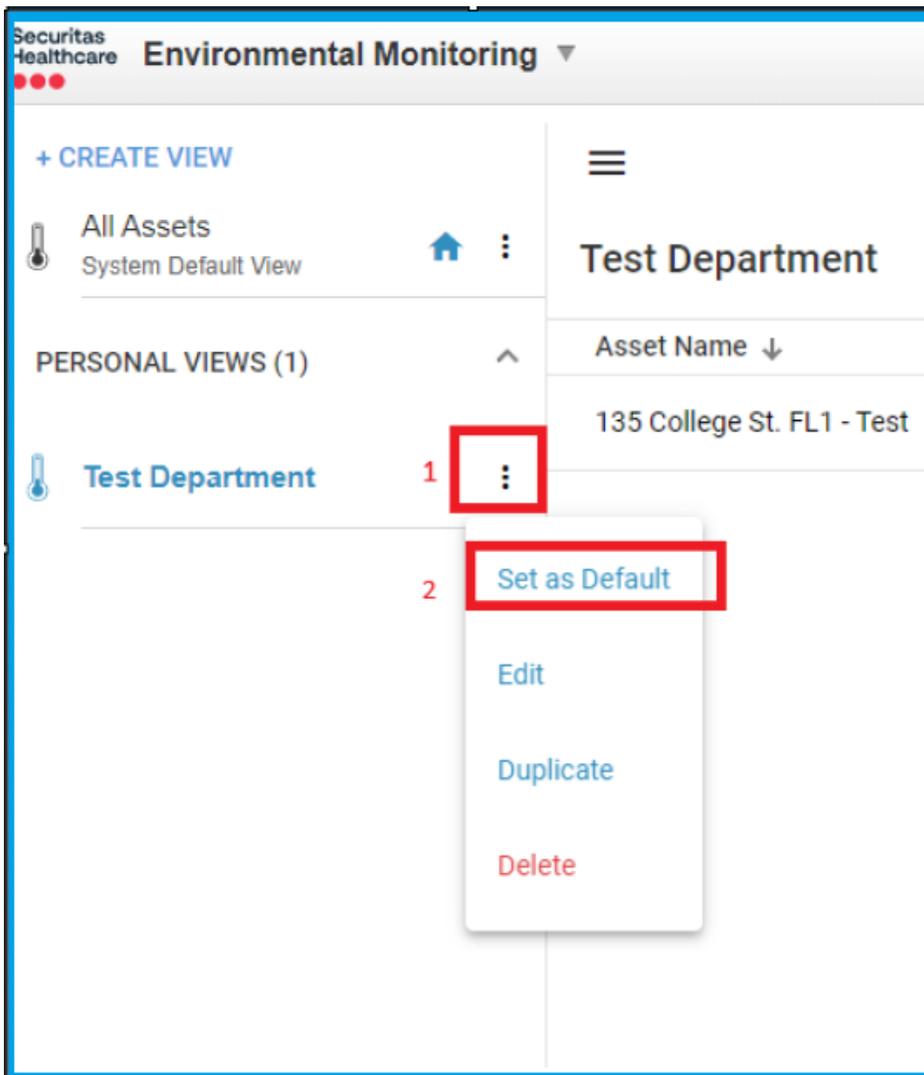

How to set up change preferences for Environmental Monitoring in MobileView Application:

To change the Time Zone, Date Format, or Temperature Units. Click the drop-down menu with the gear symbol in the upper right side of the screen. Then select My Settings. From the My Settings Menu click on the Locale tab. Select the desired preferences settings from the drop-down menu. Click OK to save settings. You must log out and back in for changes to take effect.

